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s

63

- PDQeX *forensic*GEM Sperm is a method that lyses cells and removes nucleoproteins from the DNA.
- There is no concentration step in the procedure and so the concentration of the extract is dependent on: 1) The quality of the sample; 2) In the case of swabs, the type of swab and the volume of water used to wash the swab; 3) The extraction volume (which in some cases can be scaled).
- DNA extracted using the PDQeX *forensic*GEM Sperm kit is suitable for High-Throughput Sequencing (HTS) and many types of genotyping including SNP and STR analysis as well as quantitative, multiplex and end-point PCR.
- DNA extracted using PDQeX *forensic*GEM Sperm can be quantified using a qPCR or by using fluorescent dyes like Pico Green, iQuant, Qubit assays or the like. With mixed samples, a Y-specific qPCR is recommended. Nanodrop is incompatible with PDQeX reagents.
- As with any preparative method for nucleic acid extraction, best results are obtained when samples are handled at 4°C, or on ice, before and after extraction.
- For long term storage of the extracted DNA, add TE buffer to 1x (10 mM Tris, pH 7.5, 1 mM EDTA) and store at -20°C.

Stains and Swabs

The processing of the sample will vary dependent on sample type. For liquid samples, try to keep the volume of the liquid below 10 µl. With cotton swabs, add 1/4 of the swab directly to the PDQeX cartridge. Stained fabric can be swabbed or small portions added directly to the tube.

Preparing the extraction mixture

For each extraction, make up:

10 µl 10x **ORANGE+** Buffer

2 µl *forensicGEM*

10 µl *Acrosolv*

DNA-free water - Add to a final volume of 100 µl

(Note - this will vary depending on whether you add any liquid with your sample).

PDQeX extraction

1. Add the extraction mix with sample into each PDQeX cartridge (100 µl)
2. Put the cap on the PDQeX cartridge by completely inserting the tapered column into the cartridge.
3. Load into the collection drawer either:
 - 24 well plate
 - 8 strip tubes
 - Individual tubes
4. Put the drawer in place.
5. Insert the PDQeX cartridges into the heating block.
6. Cover the cartridges with the hinged flap and close the sliding door.

MAKE SURE THE PDQeX CARTRIDGES CORRESPOND WITH A COLLECTION TUBE OR WELL- OTHERWISE YOU WILL LOSE YOUR SAMPLE

7. Select the 'Sperm' program for the extraction:

52°C 5 mins

75°C 3 mins

95°C 2 mins

- Times may be adjusted by internal laboratory optimisation.
- Changes to the default temperatures are not recommended.